

## Performa® DTR V3 96-Well Short Plate Kits

PRODUCT	CATALOG #	PURIFICATIONS
Performa DTR V3 96-Well Short Plate Kit (2 Plates)	63887	192
Performa DTR V3 96-Well Short Plate Kit (10 Plates)	89939	960

### Description

Performa DTR (**D**ye **T**erminator **R**emoval) 96-Well Plates are gel filtration plates that consist of 440 µL volume columns in a standardized array. This plate provides optimal performance for removal of unincorporated BigDye® v1.1, v3.0 and v3.1 and other dye terminators, dNTPs, salts, and other low molecular weight materials from sequencing reactions. These columns also remove DNA primers and fragments up to 20 bases, buffers, and nucleotides labeled with biotin, isotopes, and other assorted markers.

The columns are pre-packed with a fully hydrated matrix to afford optimal handling and performance characteristics. To minimize the potential for interference with sequencing applications, no preservatives, salts, or buffers are used in the preparation of these columns. Both ends of the Performa DTR 96-Well Plates are sealed to prevent drying.

The sample can be spun directly into the EdgeBio 96-Well Semi-Skirted Capillary Plates, PN 13506 (25 plates) or 90904 (100 plates), or ABI PRISM MicroAmp® Optical 96 Well Reaction Plate, thereby saving a transfer step.

COMPONENT	63887	89939
Performa DTR V3 96-Well Plate	2 plates (2x PN 4050203)	10 plates (10x PN 4050203)
Universal Plate Lids	2 lids (PN 4050094)	10 lids (2x 4050095)
96-Well Waste Plates	2 plates (PN 4050096)	10 plates (2x PN 4050097)

COMPONENT	63887	89939
96-Well Semi-Skirted Capillary Plates	2 plates (PN 4050206)	10 plates (2x PN 4050205)

### Equipment and Materials Required

1. Variable speed centrifuge (benchtop or floor model)
2. Rotor and microplate carriers for above

### Storage Condition

Store at +4°C. Do not freeze. See product label for expiration date. Alternatively, Performa DTR plates can be stored continuously at ambient temperature (defined as 20–24°C) unopened in the original packaging for up to 4 weeks from the beginning of ambient storage (the date of shipment).

### Quality Control

Field-tested for sequence quality and sequencing accuracy on capillary sequencers.

### Recommended Protocol for 5 µL–15 µL Sequencing Reaction Volumes

1. Bring reaction volume to at least 10 µL with distilled water before adding to the Performa DTR 96-Well Plate.
2. Remove the bottom and top adhesive tapes from the Performa 96-Well Plate. Cover with lid.
  - Note: Remove the bottom adhesive tape first.
  - Ensure that the plate remains horizontal to avoid losing any gel.
3. Stack the Performa 96-Well Plate on top of a 96-well waste plate. Place assembly on a cushioned centrifuge carrier.
4. Centrifuge for 3 minutes at 850 x g.<sup>1</sup> Discard eluate.
  - See “Additional Notes” for determination of RPM from RCF or visit our website at [www.edgebio.com](http://www.edgebio.com) and click on Support.
5. Transfer the reaction samples in a volume of 10–15 µL to the center of each well in the Performa 96-Well Plate. Pipet slowly. Do not touch the sides of the wells. Cover with lid.
6. Stack the Performa 96-Well Plate on top of a 96-Well Semi-Skirted Capillary Plate. Place the assembly on cushioned centrifuge carrier.

7. Centrifuge for 5 minutes at 850 x *g*. Retain eluate.
  - The eluate in the Semi-Skirted Capillary Plate contains purified sample and can be loaded directly into the DNA sequencing instrument.
  - Note: Consult the instrument manufacturer's recommendation for sample handling.

## Additional Notes

1. Conversion of RCF to RPM Calculation:

An accurate determination of the centrifugation speed is very important. The relative centrifugal force (RCF) specified in the protocol is converted to revolutions per minute (RPM) using the following formula:

$$RCF = 1.12r \left( \frac{RPM}{1000} \right)^2$$

The radius, *r*, is equal to the distance in millimeters between the axis of rotation and the bottom of the gel bed when the plate is placed in the plate carrier in the centrifuge bucket.

After measuring the radius for the specific centrifuge and accessories to be used, the proper RPM setting is calculated as follows:

$$RPM = 1000 \sqrt{\frac{RCF}{1.12r}}$$

To achieve RCF = 850 x *g*:

$$RPM = 27,549 \sqrt{\frac{1}{r}}$$

Visit the EdgeBio YouTube channel for an **RCF to RPM Conversion tutorial**.

P/N 02188 v.1

©2015 Edge BioSystems, Inc. All rights reserved. Performa® is a registered trademark of Edge BioSystems. ABI PRISM® and BigDye® are registered trademarks of Applied Biosystems or its subsidiaries in the U.S. and certain other countries.

edgebio.com

For research use only. Not for use in diagnostic procedures.



T 800-326-2685  
301-990-2685  
F 301-990-0881  
201 Perry Parkway, Suite 5  
Gaithersburg, MD 20877